



- 1 The electrode input of the PREAMP is designed for "Combination" electrodes. If separate "mono" electrodes (pH Glass and Reference) are required, specify and use BJC adapter P/N EO-1003.
- 2 When automatic temperature compensation (A.T.C.) is not required, place a 301Ω (1%) fixed resistor across PREAMP J-BOX OUTPUT terminals 1 and 2 in lieu of the A.T.C. element.
- 3 Proceed to install wiring from preamp J-Box output terminal strip to the GREAT LAKES Model 690 input terminal strip . Recommend use of multi-conductor cable as provided by GREAT LAKES, Belden #8724 or equivalent.

PARTS LIST	
DESCRIPTION	P/N
PREAMP	500-P543
JUNCTION BOX	500-P101

<b>Broadley-James Corporation</b> 1714 South Lyon Street, Santa Ana, CA 92705 Tel: (714) 547-8061 Fax: (714) 547-8988			
<b>TITLE: STANDARD PREAMP INTERFACE TO GREAT LAKES INSTRUMENTS MODEL 690</b>			
DRAWN BY: R. RAMUS	APPROVED BY:	REVISION:	DRAWING NUMBER
DATE: 10/29/91	DATE:	DATE:	<b>D-50086</b>